

N7

Low Toxicology, pH Neutral Detergent Solution



Use Instructions

InSpec N7 is designed for cleaning hard surfaces by spraying, wiping and mopping. Hold spray approximately 15cm to 20cm from the surface area. 5L versions, pour into an appropriate container for mopping. 100ml concentrate added to 2.5L of water is required for mopping. Apply to surface to ensure complete coverage.

Technical Data

N7 Ready to Use (RTU) is a 4% neutral detergent, pH7. The neutral detergent has surfactant and sequestering agents designed to assist the removal of process residue and light soiling in a cleanroom environment.

Storage

Store upright in original closed containers, away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is available in the Safety Data Sheet (SDS).

Safe Handling

Always wear gloves and goggles or face protection. Always read the label and SDS before use.

Material Compatibility

Application of solutions, when used as directed, will not affect materials normally encountered. Compatibility information in the Technical File.

Benefits

- Designed for cleaning floors, walls, ceiling and other hard surfaces for the removal of process derived residue and light soiling.
- Ideal for critical surfaces with soiling near to product manufacturing
- Very low foaming, low residue detergent with an excellent low toxicology profile
- Suitable for manual cleaning and available in convenient ready to use and concentrate formats.
- Low concentration
- 3 months validated 'in-use' shelf life
- All formats manufactured with WFI quality water
- Multiple bags for cleanroom transfer
- Manufactured to GMP

InSpec™ N7 Product

Please use the AMD Code when ordering

Product	Pack Size	Case Size	Redditch Code	AMD Code
N7 1L Manual Detergent	1L Trigger Spray	6 x 1L	N7WFI1CC-1LS	D06C10D
N7 Concentrate	100ml Screw Cap	50 x 100ml	N7WFI1-100MLS	D50C01D
N7 Session Wipes	15 piece	25 x 15pc	N7WFI1PCH15S	W25C15D
N7 Wipes	50 piece	16 x 50pc	N7WFI1PCH50S	W16C50D

AMD are the official UK distributor for Redditch Medical's range to pharmacies and compounders

Technical User Information

InSpec N7 is available in sterile RTU pre-diluted formats and as a concentrate. InSpec N7 is manufactured to GMP in an ISO 6 cleanroom. The product is filled and capped in an ISO class 5 environment.

RTU solutions are diluted with WFI quality water.

The solution is filtered through a 0.2 micron filter and gamma irradiated to a validated process to give a sterility assurance level SAL of 10⁻⁶.

InSpec N7 is supplied with Certificates of Conformity, Analysis, Irradiation and Sterility.

For screw cap bottles use the entire contents after opening.