

+ InSpec™ IPA

Sterile Isopropyl Alcohol 70% v/v with WFI Quality Water. Low residue, rapid action, fast drying, alcohol disinfectant.



+ INFORMATION

Microbiological Minimum Efficacy Testing Contact

EN 1276 Bacteria	5 minutes	EN 13697 Bacteria	5 minutes
EN 1650 Fungi	15 minutes	EN 13697 Fungi	15 minutes
EN 13704	1 minute (Bacillus Subtilis)		

+ ORDERING INFORMATION

InSpec™ IPA Products

Product	Pack Size	Case Size	Redditch Code	AMD Code
InSpec™ IPA 500ml Premium	500ml Trigger Spray	8 x 500ml	IPAWF11-500MLS	D08C50P
InSpec™ IPA 1L Premium	1L Trigger Spray	6 x 1L	IPAWF11-LS	D06C10P
InSpec™ IPA 5L Screw Cap	5 Litre Screw Cap	2 x 5L	IPAWF11-5LS	D02C00P

Please use the AMD Code when ordering.

AMD are the official UK distributor for Redditch Medical's range to pharmacies and compounders.

+ PRODUCT INFORMATION

Use Instructions

InSpec™ IPA is designed for spraying, wiping and mopping applications. Hold spray approximately 15cm to 20cm from area to be treated. 5L versions, pour into an appropriate container for mopping. Apply to surface to ensure complete coverage. For screw cap bottles use the entire contents after opening.

Technical Data

Appearance: Clear liquid free from particles. Specific Gravity (20°C): 0.872 to 0.883. Alcohol proportion: 70% +/- 2% v/v.

Storage

Store upright in original closed containers away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is available in the Safety Data Sheet (SDS).

Safe Handling

Always wear gloves and goggles or face protection. Always read the label and SDS before use.

Material Compatibility

Application of solutions, when used as directed, will not affect materials normally encountered in the cleanroom. See compatibility information in the Technical File.

Benefits

- Low residue
- Rapid action
- Bactericidal and Fungicidal
- 3 months validated 'in-use' shelf life
- Contents protected using 'Bag-in-bottle' presentation with Premium range
- Contents protected using filtered trigger presentation with standard range
- 1L and 5L formats manufactured with WFI quality water
- Multiple bags for cleanroom transfer
- Ideal surface and aseptic transfer disinfectants
- Manufactured to GMP

Technical User Information

InSpec™ IPA is a sterile ready to use disinfectant. This is a 70% v/v alcohol solution in WFI quality water.

InSpec™ IPA is manufactured to GMP in an ISO 6 cleanroom. The product is filled and capped in an ISO class 5 environment.

The solution is filtered through a 0.2 micron filter and gamma irradiated to a validated process (at 25-45kGy) to give a sterility assurance level SAL of 10⁻⁶.

InSpec™ IPA is supplied with Certificates of Conformity, Analysis, Irradiation Sterility and Endotoxin.

All formats of InSpec™ IPA have a stabilised 24 month shelf life.

Redditch Medical

Discovery 2, NETPark, 2 William Armstrong Way, Sedgefield, County Durham, TS21 3FD
UK Tel: +44 (0) 1332 755622 | International Tel: +44 (0) 1527 830942
sales@redditchmedical.com | www.redditchmedical.com