

## + InSpec® IPA

Sterile Isopropyl Alcohol, Sterile Wipes, 70% v/v with WFI Quality Water.

Low residue, rapid action, fast drying, alcohol disinfectant.



## + INFORMATION

### Microbiological Minimum Efficacy Testing Contact

EN 1276 Bacteria	5 minutes	EN 13697 Bacteria	5 minutes
EN 1650 Fungi	15 minutes	EN 13697 Fungi	15 minutes

## + ORDERING INFORMATION

### InSpec® IPA Products

Product	Pack Size	Case Size	Redditch Code	AMD Code
InSpec® IPA 500ml Premium	500ml Trigger Spray	8 x 500ml	IPAWF11-500MLS	D08C50P
InSpec® IPA 1L Premium	1L Trigger Spray	6 x 1L	IPAWF11-LS	D06C10P
InSpec® IPA 5L Screw Cap	5 Litre Screw Cap	2 x 5L	IPAWF11-5LS	D02C00P
InSpec® IPA Pouch 15pc	15 Wipes	25 per case	IPAWFIPCH15S	W25C15P
InSpec® IPA Pouch 50pc	50 Wipes	16 per case	IPAWFIPCH50S	W16C50P
InSpec® IPA Standard Tub	200 Wipes	12 per case	IPAWF112TB200S	W12C20P

Please use the AMD Code when ordering.

AMD are the official UK distributor for Redditch Medical's range of products to the NHS, pharmacies and compounders.

## + PRODUCT INFORMATION

### Use Instructions

InSpec® IPA is designed for spraying, wiping and mopping applications. Hold spray approximately 15cm to 20cm from area to be treated. 5L versions, pour into an appropriate container for mopping. Apply to surface to ensure complete coverage. For screw cap bottles use the entire contents after opening.

### Technical Data

Appearance: Clear liquid free from particles. Specific Gravity (20°C): 0.872 to 0.883. Alcohol proportion: 70% +/- 2% v/v.

### Storage

Store upright in original closed containers away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is available in the Safety Data Sheet (SDS).

### Safe Handling

Always wear gloves and goggles or face protection. Always read the label and SDS before use.

### Material Compatibility

Application of solutions, when used as directed, will not affect materials normally encountered in the cleanroom. See compatibility information in the Technical File.

### Benefits

- Low residue
- Rapid action
- Bactericidal and Fungicidal
- Contents protected using 'Bag-in-bottle' presentation with Premium range
- Contents protected using filtered trigger presentation with standard range (spray bottle)
- 1L and 5L formats manufactured with WFI quality water
- Multiple bags for cleanroom transfer
- Ideal surface and aseptic transfer disinfectants
- Manufactured to GMP
- Active Substance compliant with Article 95 of the Biocidal Products Regulation (BPR)

### Technical User Information

InSpec® IPA is a sterile ready to use disinfectant. This is a 70% v/v alcohol solution in WFI quality water.

InSpec® IPA is manufactured to GMP and is filled and capped in an ISO class 5 environment.

The solution is filtered through a 0.2 micron filter and gamma irradiated to a validated process (20-50kGy for 5L products, 25-45kGy for all other sizes) to give a sterility assurance level SAL of 10<sup>-6</sup>.

### Aseptic Medical Devices

Newmarket Drive, Derby DE24 8SW | +44 (0) 1332 755622  
sales@asepticmedical.com | www.asepticmedical.com

Aseptic Medical Devices is a trading name of Riverside Medical Packaging Company Limited