

+ InSpec™ N10

Manual Detergent Pre-diluted. An alkaline detergent and hard surface cleaner for use in the biopharmaceutical and cleanroom industries.



+ ORDERING INFORMATION

InSpec™ N10 Products

Product	Pack Size	Case Size	Redditch Code	AMD Code
InSpec™ N10 100ml Detergent Sterile Concentrate	100ml screw cap	50 x 100ml	N10WF11-100MLS	D50C01N
InSpec™ N10 1L Manual Detergent Sterile RTU	1L trigger spray	6 x 1L	N10WF11CC-1LS	D06C10N

Please use the AMD Code when ordering.

AMD are the official UK distributor for Redditch Medical's range to pharmacies and compounders.

+ PRODUCT INFORMATION

Use Instructions

InSpec™ N10 is designed for cleaning hard surfaces by spraying, wiping and mopping. Hold spray approximately 15cm to 20cm from the surface area. 5L versions, pour into an appropriate container for mopping. 100ml concentrate added to 2.5L of water is required for mopping. Apply to surface to ensure complete coverage.

Technical Data

InSpec™ N10 Ready to Use (RTU) is a 4% alkaline detergent, pH12. The concentrate presentation has a pH of >13. The alkaline detergent has surfactant and sequestering agents designed to assist the removal of process residue and general soiling.

Storage

Store upright in original closed containers, away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is available in the Safety Data Sheet (SDS).

Safe Handling

Always wear gloves and goggles or face protection. Always read the label and SDS before use.

Material Compatibility

Application of solutions, when used as directed, will not affect materials normally encountered in the cleanroom. See compatibility information in the Technical File.

Benefits

- Designed for cleaning floors, walls, ceiling and other hard surfaces for the removal of general process derived residue and soiling
- The Dycem recommended cleaner of Dycem flooring
WFI quality water
- Multiple bags for cleanroom transfer
- Manufactured to GMP
- Rapid action alkaline detergent. Suitable for manual cleaning. Available in convenient ready to use formats
- 3 months validated 'in-use' shelf life
- 1L and 5L formats manufactured with WFI quality water

Technical User Information

InSpec™ N10 is available as a sterile and non-sterile detergent in pre-diluted and concentrated formats.

InSpec™ N10 is manufactured to GMP in an ISO 6 cleanroom. The product is filled and capped in an ISO class 5 environment.

RTU solutions are diluted with WFI quality water.

The solution is filtered through a 0.2 micron filter and gamma irradiated to a validated process (at 25-45kGy) to give a sterility assurance level SAL of 10⁻⁶.

InSpec™ N10 is supplied with Certificates of Conformity, Analysis, Irradiation and Sterility.

For screw cap bottles use the entire contents after opening.

Redditch Medical

Discovery 2, NETPark, 2 William Armstrong Way, Sedgefield, County Durham, TS21 3FD
UK Tel: +44 (0) 1332 755622 | International Tel: +44 (0) 1527 830942
sales@redditchmedical.com | www.redditchmedical.com