

+ InSpec™ QT+

Sterile Quaternary Ammonium Compound, Disinfectant Cleaner. Exceptional Quat disinfectant with proven efficacy against a wide range of Bacteria, Yeast and Fungi.



+ INFORMATION

Microbiological Minimum Efficacy Testing Contact

EN 1276 Bacteria	5 minutes	EN 13697 Bacteria	5 minutes
EN 1650 Fungi	15 minutes	EN 13697 Fungi	15 minutes

+ ORDERING INFORMATION

InSpec™ QT+ Products

Product	Pack Size	Case Size	Redditch Code	AMD Code
InSpec QT+ Sterile Concentrate 102ml	1	50 x 102ml	QTPLCNC-102	D50C1QT
InSpec QT+ Sterile 900ml Trigger RTU	1	6 x 900ml	QTPLWFI31-900MLS	D06C9QT
InSpec QT+ Sterile 5 L RTU	1	2 x 5L	QTPLWFI31-5LS	D02C0QT

Please use the AMD Code when ordering.

AMD are the official UK distributor for Redditch Medical's range to pharmacies and compounders.

+ PRODUCT INFORMATION

Use Instructions

InSpec QT+ is designed for spraying, wiping and mopping applications. Hold spray approximately 15cm to 20cm from area to be treated. 5L versions, pour into an appropriate container for mopping. 102ml concentrate added to 5.0L of water is required for mopping. Apply to surface to ensure complete coverage.

Technical Data

Formulated Quaternary Ammonium Compound based disinfectant/cleaner. Active ingredient 0.12% RTU (ready to use). RTU pH = 10.5 to 11.5. Concentrate pH = 12.5 to 13.5.

Storage

Store upright in original closed containers, away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is available in the Safety Data Sheet (SDS).

Safe Handling

Always wear gloves and goggles or face protection. Always read the label and SDS before use.

Material Compatibility

Application of solutions, when used as directed, will not affect materials normally encountered in the cleanroom. See compatibility information in the Technical File.

Benefits

- An exceptional non-oxidising biocide
- Bactericidal and Fungicidal
- Low residue
- Contents protected using 'bag-in-bottle' technology. Provides a validated 3 month 'in-use' shelf life.
- 1L and 5L formats manufactured with WFI quality water
- Double bagged to enable rapid cleanroom transfer
- Ideal rotational partner for InSpec™ AN, InSpec™ HA and OX products
- Manufactured in accordance with GMP

Technical User Information

InSpec QT+ is a sterile disinfectant in pre-diluted and concentrated formats.

InSpec QT+ is manufactured in accordance with GMP in an ISO 6 cleanroom. The product is filled and capped in an ISO class 5 environment.

The solution is filtered through a 0.2 micron filter and gamma irradiated via a validated process (at 25-45kGy) providing a sterility assurance level SAL of 10⁻⁶.

Trigger sprays supplied with 900ml solution in a 1L container.

InSpec QT+ is supplied with certificates of Conformity, Analysis, Irradiation and Sterility. All formats of InSpec QT+ have a stabilised 24 month shelf life.

For screw cap bottles use the entire contents after opening.

Redditch Medical

Discovery 2, NETPark, 2 William Armstrong Way, Sedgefield, County Durham, TS21 3FD
UK Tel: +44 (0) 1332 755622 | International Tel: +44 (0) 1527 830942
sales@redditchmedical.com | www.redditchmedical.com