

## + InSpec® QT

Sterile Quaternary Ammonium Compound, Disinfectant Cleaner. Exceptional Quat disinfectant showing anti-microbial properties. Proven efficacy against a wide range of bacteria, yeast and mould.

## + INFORMATION

### Microbiological Minimum Efficacy Testing Contact

EN 1276 Bacteria	5 minutes	EN 13697 Bacteria	5 minutes
EN 1650 Fungi	15 minutes	EN 13697 Fungi	15 minutes

## + ORDERING INFORMATION

### InSpec® QT Products

Product	Pack Size	Case Size	Redditch Code	AMD Code
InSpec® QT 900ml Trigger	900ml Trigger Spray	6 x 900ml	QTWFI30-1LS	D06C90Q
InSpec® QT 5L Screw Cap	5 Litre Screw Cap	2 x 5L	QTWF130-5LS	D02C00Q
InSpec® QT Concentrate	100ml Screw Cap	50 x 100ml	QTCNC30-100	D50C01Q
InSpec® QT Burstable Pouch	15 wipes/pouch	20 per case	QTPLWFI31BP15S	W10C2QT

Please use the AMD Code when ordering.

AMD are the official UK distributor for Redditch Medical's range of products to the NHS, pharmacies and compounders.



## + PRODUCT INFORMATION

### Use Instructions

InSpec® QT is designed for spraying, wiping and mopping applications. For spraying hold approximately 15 to 20cm from area to be treated. For screw cap versions, pour into an appropriate container for mopping and add 100ml of concentrate per 2.4L of water. Apply to surface to ensure complete coverage.

### Technical Data

Aldehyde-free disinfectant cleaner. RTU pH 11.5 to 12.5, concentrate pH 12.5.

### Storage

Store upright in original closed containers away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is available in the Safety Data Sheet (SDS).

### Safe Handling

Always wear gloves and goggles or face protection. Always read the label and SDS before use.

### Material Compatibility

Application of solutions, when used as directed, will not affect materials normally encountered in the cleanroom. See compatibility information in the Technical File.

### Benefits

- An exceptional non-oxidising biocide
- Bactericidal and Fungicidal
- Cleans and disinfects in the presence of dirt, blood and proteins.
- Contents protected using 'bag-in-bottle' presentation with premium range (spray bottle)
- 1L and 5L formats manufactured with WFI quality water
- Multiple bags for cleanroom transfer
- Ideal rotational partner for InSpec® AN, InSpec® HA and InSpec® OX
- Manufactured to GMP
- Active Substance compliant with Article 95 of the Biocidal Products Regulation (BPR)

### Technical User Information

InSpec® QT is a sterile disinfectant in pre-diluted and concentrated formats.

InSpec® QT is manufactured to GMP and is filled and capped in an ISO class 5 environment.

The solution is filtered through a 0.2 micron filter and gamma irradiated to a validated process (at 20-50kGy for 5L products and 25-45kGy for all other formats) to give a sterility assurance level SAL of 10<sup>-6</sup>.

Trigger sprays supplied with 900ml solution in a 1L container.

For screw cap bottles use the entire contents after opening.

### Aseptic Medical Devices

Newmarket Drive, Derby DE24 8SW | +44 (0) 1332 755622  
sales@asepticmedical.com | www.asepticmedical.com

Aseptic Medical Devices is a trading name of Riverside Medical Packaging Company Limited