

+ InSpec™ QT

Sterile Quaternary Ammonium Compound, Disinfectant Cleaner. Exceptional Quat disinfectant showing anti-microbial properties. Proven efficacy against a wide range of bacteria, yeast and mould.



+ INFORMATION

Microbiological Minimum Efficacy Testing Contact

EN 1276 Bacteria	5 minutes	EN 13697 Bacteria	5 minutes
EN 1650 Fungi	15 minutes	EN 13697 Fungi	15 minutes

+ ORDERING INFORMATION

InSpec™ QT Products

Product	Pack Size	Case Size	Redditch Code	AMD Code
InSpec™ QT 900ml Trigger	900ml Trigger Spray	6 x 900ml	QTFI30-1LS	D06C90Q
InSpec™ QT 5L Screw Cap	5 Litre Screw Cap	2 x 5L	QTFI30-5LS	D02C00Q
InSpec™ QT Concentrate	100ml Screw Cap	50 x 100ml	QTCNC30-100	D50C01Q

Please use the AMD Code when ordering.

AMD are the official UK distributor for Redditch Medical's range to pharmacies and compounders.

+ PRODUCT INFORMATION

Use Instructions

InSpec™ QT is designed for spraying, wiping and mopping applications. Hold spray approximately 15cm to 20cm from area to be treated. 5L versions, pour into an appropriate container for mopping. 100ml concentrate added to 2.5L of water is required for mopping. Apply to surface to ensure complete coverage.

Technical Data

Aldehyde-free disinfectant cleaner.
4% RTU Solution pH 11.8-12.2
Concentrate pH 12.9.

Storage

Store upright in original closed containers away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is available in the Safety Data Sheet (SDS).

Safe Handling

Always wear gloves and goggles or face protection. Always read the label and SDS before use.

Material Compatibility

Application of solutions, when used as directed, will not affect materials normally encountered in the cleanroom. See compatibility information in the Technical File.

Benefits

- An exceptional non-oxidising biocide
- Bactericidal and Fungicidal
- Clean and disinfects in the presence of dirt, blood and proteins.
- 3 months validated 'in-use' shelf life
- Contents protected using 'bag-in-bottle' presentation with premium range
- 1L and 5L formats manufactured with WFI quality water
- Multiple bags for cleanroom transfer
- Ideal rotational partner for InSpec™ AN, InSpec™ HA and InSpec™ OX
- Manufactured to GMP

Technical User Information

InSpec™ QT is a sterile disinfectant in pre-diluted and concentrated formats. InSpec™ QT is manufactured to GMP in an ISO 6 cleanroom.

The product is filled and capped in an ISO class 5 environment.

The solution is filtered through a 0.2 micron filter and gamma irradiated to a validated process (at 25-45kGy) to give a sterility assurance level SAL of 10⁻⁶.

Trigger sprays supplied with 900ml solution in a 1L container.

InSpec™ QT is supplied with Certificates of Conformity, Analysis, Irradiation Sterility and Endotoxin.

All formats of InSpec™ QT have a stabilised 24 month shelf life.

For screw cap bottles use the entire contents after opening.

Redditch Medical

Discovery 2, NETPark, 2 William Armstrong Way, Sedgefield, County Durham, TS21 3FD
UK Tel: +44 (0) 1332 755622 | International Tel: +44 (0) 1527 830942
sales@redditchmedical.com | www.redditchmedical.com